Manual de instrucciones User manual

Ultrasonidos Estmon
Estmon Ultrasonic Cleaners

Ref. LIM0E0220 - 2,2 L

Ref. LIM0E0300 - 3,0 L

Ref. LIM0E0450 - 4,5 L

Ref. LIM0E0600 - 6,0 L

Ref. LIM0E1500 - 15 L

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1. Introduction

*Congratulations on the purchase of your new **B Estmon** ULTRASONIC CLEANER ELECTRONIC SERIES. You have just entered the world of smart cleaning!*

Following these instructions will allow you to use this ultrasonic cleaning equipment, taking full advantage of its advanced technology.

It is a device manufactured with high quality electronic components, it is equipped with a fully digital multifunction LCD control panel. This allows us to visualize and control by means of the touch keyboard all the necessary parameters to obtain an excellent ultrasonic cleaning.

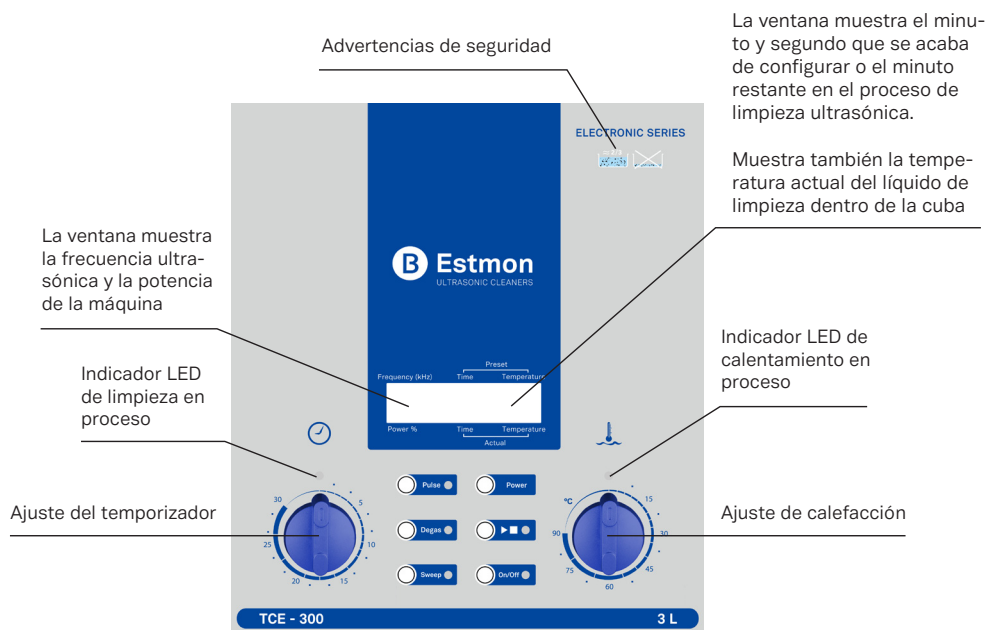
Among its most advanced functions we can highlight the multidirection sweep function and the power increase function up to 120%, increasing efficiency and equipment performance.

Fully manufactured under the most demanding European standards, with CE certificate.

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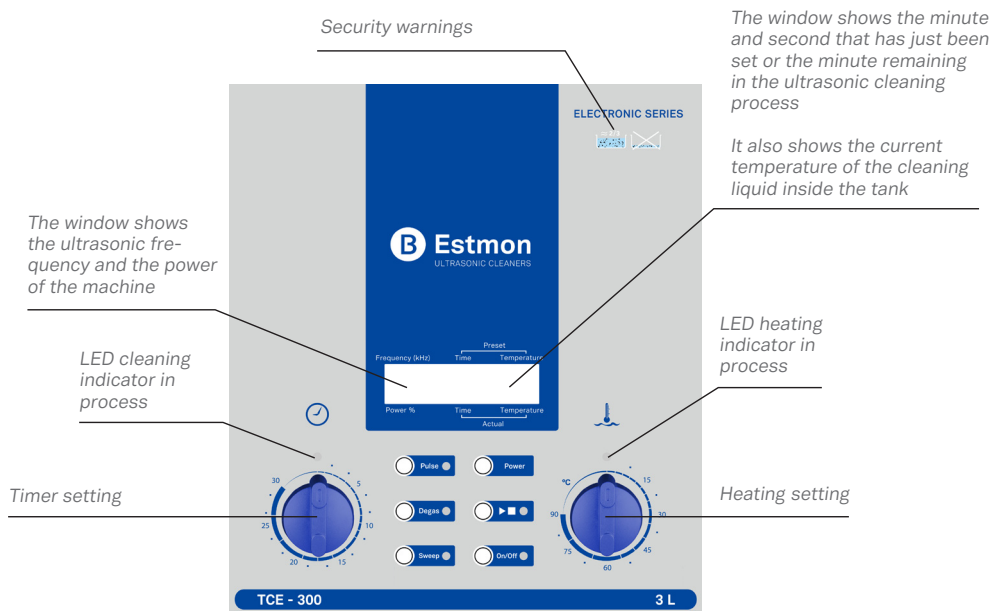
2. Use Guide

2.1 Machine controls

Legend of buttons and symbols on the screen

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2.2 Start up

- *Before turning on the power, place the unit on a stable and flat work surface. Add water to the tank, suitable to completely submerge the items to be cleaned. It is recommended to use a stainless steel basket.*

*Never use the unit without water in the cleaning tank.
Always fill the unit above the 2/3 marker line indicated
inside the tank.*

Otherwise, DAMAGE MAY OCCUR.

- *Connect the power cord to the back of the unit and make sure the wall supply is ON. All buttons are touch sensitive, just lightly touch with your finger. Touch the ON / OFF button on the front panel to turn on the machine.*

In normal operating state the power is 100% and the ultrasonic frequency is 37 kHz.

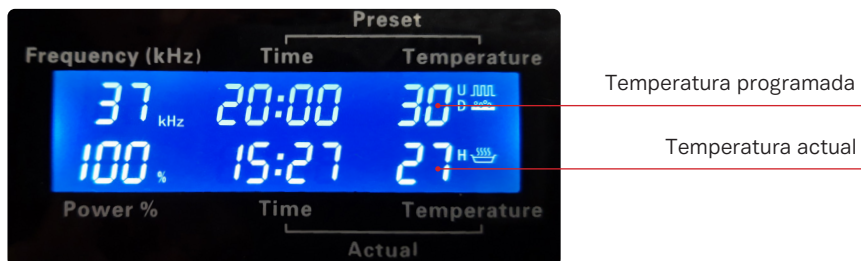
2.3 Temperature setting

- *When turned on, the LCD display module shows the actual water temperature (Actual temperature). Turn the heating knob clockwise to freely set the required temperature, this can be adjusted by 5 degree steps:*

30 - 35 - 40 - 45 - 50 - 55 - 60 - 65 - 70 - 75 - 80 °C

Once the temperature is set, the green LED located on the top of the control will illuminate.

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3 - 5 - 8 - 10 - 13 - 15 - 18 - 20 - 23 - 25 - 30 min.

The optimum working temperature is 60 °C.

When the liquid in the tank reaches the selected temperature, the heating will stop automatically and the symbol H: ■ appears. If the temperature is below the set temperature, the heating will also start automatically, appearing  symbol.

- The heating function is not restricted by the ultrasonic cleaning time, although it can work for a long time (however, it is necessary to ensure that there is enough water with respect to the level marked in the tank).

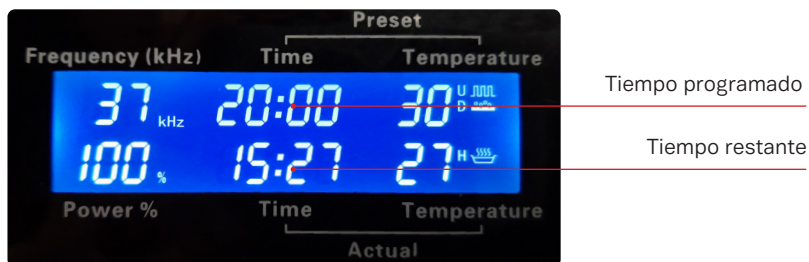
2.4 Timer setting

- With the machine running, turn the timer knob clockwise to set the necessary parameters. The time can be regulated by different periods of time, comprising:

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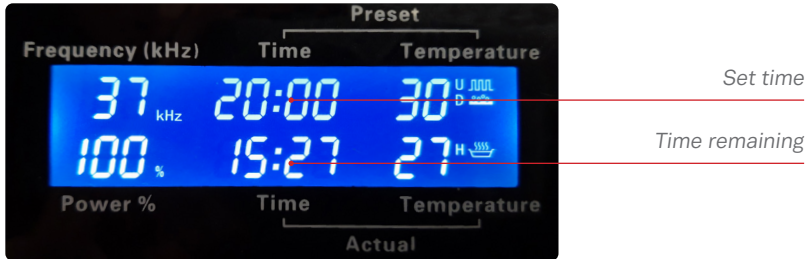
Once the time is set, the green LED located on the top of the control will light up.

The LCD module will show the pre-selected settings. At this time, the ultrasonic cleaning function will start quickly if the ►■ button is pressed. The LCD module will show the remaining time.

When the cleaning time is over, the ultrasound will stop. If we want to repeat the cleaning cycle, we only have to press the ►■ button. Cleaning will start again.

2.5 Power regulation options

- The factory default settings for the machine are programmed with 100% power. The power range varies from 40% to 70% and 100%, which can be freely chosen by pressing the Power button. Then the set value will be displayed on the LCD module.

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2.5.1 Ultrasonic power increase function

- Once the *Pulse* button is pressed, the machine power will increase to 120% and the button indicator light will turn on. If the *Pulse* button is pressed again, the power increase function will stop. The indicator light on the button will turn off and the power will resume the standard value.

2.6 Ultrasonic wave options

2.6.1 Degassing function

- With the  button pressed, once the *Degas* button is pressed, the machine will start the degassing function and the indicator light will come on, emitting an intermittent sound that will decrease as the process is completed.

Degas: D 

It is important to perform this function with a minimum period of 15 minutes, and whenever the water is changed. This will offer less resistance to ultrasonic waves. This will have a positive impact on an increase in power and performance of the machine.

Once the *Degas* button is pressed again, the degassing function will stop and the degassing button indicator light will turn off.

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2.6.2 Sweep function

- The sweep function modifies the trajectory of the ultrasonic wave creating a cross sweep that will have an impact on cleaning efficiency in areas with extreme dirt and difficult access.

This function will be activated once we have pressed the ►■ button. Next, we keep the Sweep button pressed for 2 seconds.

Sweep: S 

If the Sweep button is pressed again, the machine will stop the sweep function and the indicator light will turn off.

3. Warranty

All **B Estmon** models are guaranteed for 12 months against any manufacturing defect.

3.1 Exclusions of warranty coverage

- The improper use or manipulation of the equipment.
- The use of caustic agents or acids in the tank.
- Use of the machine with an insufficient level of water inside the tank.

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5. Notes

- *For a MORE EFFICIENT CLEANING, add the Technoflux ultrasonic cleaner 909 and 707 degreasing salts (Ref. LIM0P0200 y Ref. LIM0P0100 respectively). Other alkaline cleaning agents can also be used.*

Gabriel Benmayor S.A.
A-60512100
Bach, 2-B. Pol. Ind. Foinvasa
08110 Montcada i Reixac, Barcelona
benmayor@benmayor.com
T +34 935 724 161 / F +34 935 724 165

benmayor.com